


## User Manual – Count

- All the 4 DAQ machines will have a folder – Count in the desktop
- Count folder will have two sub folders named as – a) Stretcher\_Channel\_Count and b) Panel\_Tube\_Count
- Stretcher\_Channel\_Count reports the count for a stretcher - channel combination
- Panel\_Tube\_Count reports the count for a Panel-Tube combination
- Both the programs have similar user interface.

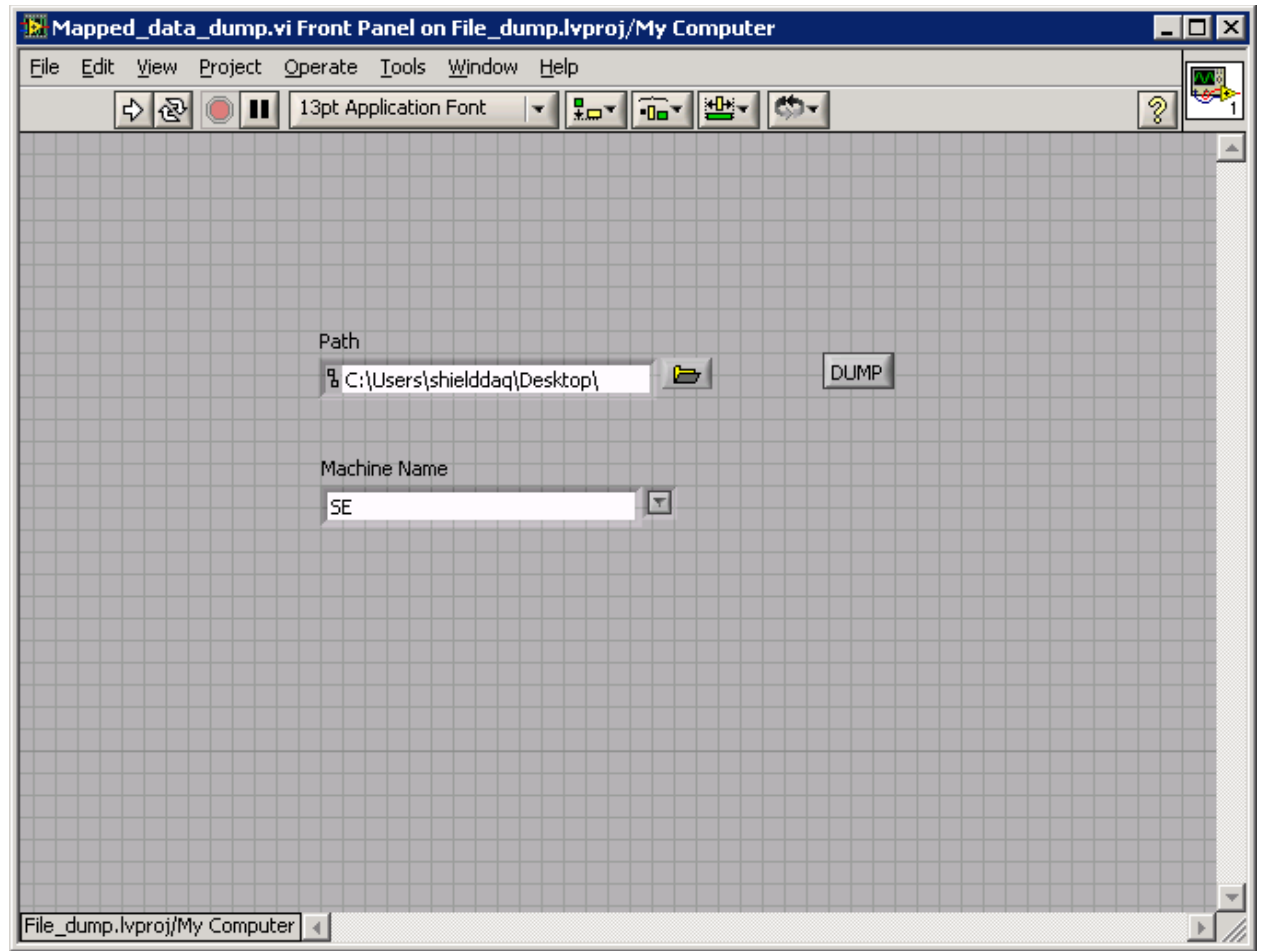

- Click on the folder icon to choose the name and path of the file to which data should be dumped.
- Choose the machine name on which you are running this program
- Click Dump button.
- In few seconds, the program stops and the data will be dumped to the file specified.

Stretcher\_Channel\_Count will have an output similar to the below image.

[illegible]

- Each column corresponds to a stretcher and each row corresponds to a channel in that stretcher.

Panel\_Tube\_Count will have an output similar to the below image.

[illegible]